

# TMB ELISA Substrate

<b>Product code:</b>	3652-F10
<b>Contents:</b>	Contains TMB, 3,3',5,5' tetramethylbenzidine, with hydrogen peroxide included.
<b>Applications:</b>	Ready-to-use substrate for ELISA assays with horseradish peroxidase.
<b>Instructions:</b>	Add the substrate after incubation with horseradish peroxidase and subsequent washing. Incubate for typically 5-15 minutes and then stop the enzymatic reaction by adding an acid, e.g. 1 M H <sub>2</sub> SO <sub>4</sub> . This will change the color from blue to yellow. Read the absorbance at 450 nm.
<b>Storage:</b>	Shipped at ambient temperature. On arrival store refrigerated at 4-8°C. Substrate should preferably be used cold.
<b>Size:</b>	120 ml



Distribuito in ITALIA da  
**Li StarFish S.r.l.**  
Via Cavour, 35  
20063 Cernusco S/N (MI)  
telefono 02-92150794  
fax 02-92157285  
info@listarfish.it  
www.listarfish.it

Note; for research use only.  
Mabtech shall not be liable for the use or handling  
of the product or for consequential, special,  
indirect or incidental damages therefrom.

